

## Auto Melting Point Apparatus

1935 & 1934

### Advantages

- Easy to Operate
- 4 Digit Seven Segment LED Display
- Programmable Plateau Temperature
- Programmable Ramp Rates
- Storage upto 100 Results
- Quick Cooling with Integrated Fan
- Audio Visual Indicators for System Status

# Auto Melting Point Apparatus



1 Year Warranty

Auto Melting Point Apparatus is specially designed to find the melting point temperature of crystalline sample held within capillary tubes. Instrument uses the latest microcontroller technology and advanced engineering techniques so as to give enhanced accuracy and reproducibility.

The system is equipped with microcontroller based digital display for displaying accurate results and soft touch membrane type keys for ease of operation.

The sample capillary to be analyzed is held in an heating block (i.e. aluminium block) inside the heating chamber. The block holding the sample capillary tube is heated and the temperature is displayed on the 4 digit seven segment red LED display.

The plateau temperature and the heating ramp rates are fully programmable using 'Plateau'/'Ramp' keys along with the 'Up', 'Down' keys. The plateau temperature allows the user to set a temperature to few degrees below the expected melting point temperature. Auto melting point apparatus will then heat to the programmed plateau temperature very rapidly. Once plateau temperature is reached the instrument will heat slowly from the plateau temperature until the melt occurs on programmed ramp rate.

The end point of the measurement cycle (i.e. final melting point temperature) is indicated with 'END' led along with the intermittent audible beep. 'Run' led indicates that the instrument is in measurement cycle

User can store upto 100 results (i.e. melting point temperatures) using the 'Store/Recall' key which are retained in the instruments memory even when it is switched off.

If by chance the sample capillary is broken inside the heating chamber, the flush the broken capillary using the 'Normal'/'Clean' rotary switch provided side of the instrument





## Auto Melting Point Apparatus

### Technical Data

<b>NUMBER OF SAMPLES</b>	1 Sample at a Time
<b>TEMPERATURE RANGE</b>	Ambient to 300 °C
<b>TEMPERATURE ACCURACY</b>	± 1 °C at 20 °C ± 2.5 °C at 300 °C
<b>TEMPERATURE SENSOR</b>	Platinum Resistance Pt100
<b>RAMP RATE</b>	Programmable at Scale of 1 – 10
<b>PLATEAU TEMPERATURE</b>	Programmable
<b>DISPLAY</b>	4 Digit Seven Segment LED Display (Model – 1935) 3 Digit Seven Segment LED Display (Model – 1934)
<b>DISPLAY RESOLUTION</b>	0.1 °C (Model – 1935) 1 °C (Model – 1934)
<b>KEYBOARD</b>	Soft Touch Membrane Keys
<b>STATUS LED'S</b>	PLATEAU, °C, RAMP, RUN, END, READY & RECALL
<b>MEMORY</b>	Storage upto 100 Results
<b>POWER</b>	AC 230 V ± 10%, 50 Hz
<b>WEIGHT</b>	9.0 Kg. (Approx.)
<b>ACCESSORIES</b>	Box of Capillaries, Operation Manual & Dust Cover

Specifications are subject to change due to continuous

### Product Range

Semi Auto Analyser, Double Beam UV/VIS Spectrophotometer, Single Beam UV/VIS Spectrophotometer, Flame Photometer, Hemoglobin Meter, Photo Colorimeter, pH Meter, Conductivity Meter, TDS Meter, DO Meter, Salinity Meter, Turbidity Meter, Nephelometer, Colony Counter, Karl Fischer Moisture Titrator, Photo Fluorometer, Dissolution Test Apparatus, Disintegration Test Apparatus, Friability Test Apparatus, Melting Point Apparatus, Telethermometer, Temperature Indicator, Water & Soil Analysis Kit, Portable pH Meter, Portable Conductivity Meter, Portable TDS Meter, Portable DO Meter, Student Microscope, Binocular microscope, Co-Axial Binocular, Medical Microscope

